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### **Stratagene standard protocol**

1. Pre-chill a 14-ml BD Falcon polypropylene round-bottom tubes on ice. Preheat NZY+ broth to  $42^\circ\text{C}$ .
2. Thaw the cells on ice. When thawed, gently mix and aliquot  $100 \mu\text{l}$  of cells into the pre-chilled tubes.
3. Add  $4 \mu\text{l}$  of the  $\beta$ -ME ( $\beta$  巯基乙醇) to the aliquot of cells.
4. Swirl the tubes gently. Incubate the cells on ice for 10 minutes, swirling gently every 2 minutes.
5. Add  $0.1$ - $50 \text{ ng}$  of the experimental DNA (or  $2 \mu\text{l}$  of a ligation mixture) to the aliquot of cells.
6. Swirl the tubes gently, then incubate the tubes on ice for 30 minutes.

7. Heat-pulse the tubes in a 42°C water bath for 30 seconds. The duration of the heat pulse is critical.
8. Incubate the tubes on ice for 2 minutes.
9. Add 0.9 ml of preheated (42°C) NZY+ broth and incubate the tubes at 37°C for 1 hour with shaking at 225-250 rpm.
10. Plate  $\leq 200$   $\mu$ l of the transformation mixture on LB agar plates containing the appropriate antibiotic (and containing IPTG and X-gal if color screening is desired).
11. Incubate the plates at 37°C overnight. If performing blue-white color screening, incubate the plates at 37°C for at least 17 hours to allow color development (color can be enhanced by subsequent incubation of the plates for 2 hours at 4°C).